**BOMBA DE VACIO /
ROTARY VANE VACUUM PUMP
Models: XZ-1B & 2XZ-1G**

Cod. 5900620

Cod. 5900621

MANUAL DE INSTRUCCIONES /

INSTRUCTION MANUAL

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or COMECTA's technical service.*
- 5) **IMPORTANT! COMECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
FAILURES
MAINTENANCE
SPARE PARTS
GUARANTEE

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The standard equipment consist of the following components:

Pieza	Piece	Código/Code	Cant. / Qty.
Bomba Vacío XZ-1B	Vacuum Pump XZ-1B	5900620	1
Bomba Vacío 2XZ-1G	Vacuum Pump 2XZ-1G	5900621	1
Manual de Instrucciones	User Manual		1

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1. Direct vacuum connection, high speed, small volume and light weight.
2. Energy saving, easy to operate.
3. Single phase motor with 220V or 110V power supply.

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4. Low noise, low vibrations, practical and easy to use.

5. Equipped with a Gas-ballast valve to prolong the life of the pump oil.

6. The pump will operate continuously in an ambient temperature of 5°C to 40°C. The inlet pressure is less than 1330 Pa.

7. Applications suitable for Laboratory and medical applications such as degassing, rotary evaporation, centrifugal evaporation, distillation, electrophoresis gel drying, vacuum oven, chemical industry, gas analysis and general vacuum applications.

8. The motor has a termic device for protecting the apparatus when it reaches to a certain temperature. It starts on when the temperature goes down.

INSTALLATION

Place the pump on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

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Do not use the apparatus to dry or for thermal treatment of products which can give off vapours capable of making an explosive mixture.

THE PUMP IS NOT SUITABLE FOR DECATING (MIXING) LIQUIDS.

A VACUUM TRAP IS HIGHLY RECOMMENDABLE IN THE CONNECTION TO THE PUMP TO AVOID THE ACCIDENTAL LEAK OF LIQUIDS.

Be sure that the voltage supply is the same as

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OPERATING METHOD

- 1) *Before starting to use the new Vacuum pump, you must add clean vacuum pump oil into the pump. Remove the oil seal lid. The oil must be added to the level marked on the window of the pump, not too much and not less than the marked line. Otherwise, the pump will be damaged or oil will be sprayed out if too much is added.*
- 2) *Please use the same vacuum pump oil when topping up, do not mix differing branded oils.*
- 3) *The pump should not be run in direct air for more than 3 minutes, (i.e. not connected to the item being evacuated)*
- 4) *The pump should not be used to evacuate gases with corrosive properties, dust, high humidity or products with combustible or explosive properties.*
- 5) *The tube diameter must not be larger than the inlet diameter of the pump. To reduce the possibility of the tube doubling and reducing airflow the connection should be as short and as direct as possible.*
- 6) *Make sure that all tubes and connectors are well sealed.*
- 7) *If the gas being pumped is health hazard, please connect the appropriate filter to the outlet. Exhaust outlet, with a hose to an external source, far from the working environment.*
- 8) *Connecting to the mains supply, check the voltage power rating as stipulated on the nameplate of the pump, i.e. 110V or 220V.*

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MAINTENANCE

1. The vacuum pump should be used in a dry and clean environment.
2. For new vacuum pumps, you are requested to change the oil after working 250 - 300 hours and to change the oil 600 - 1000 hours thereafter. To check the oil level, it must be done when the pump is cold.
More frequent oil changes will be required depending on the application.
3. To change the oil, run the pump for 30 minutes. Remove the oil drain valve and drain the pump. Flush the pump with fresh oil. Replace the drain valve. Then refill with oil to the mark in the window center. Replace the oil fill cap.
4. The vacuum degree will be down after the pump runs for a long time. Remove the oil box, clean the filter on the cover of the stator, change the vacuum oil and install the oil box back. Then check the vacuum degree. If it does not work, the vacuum pump needs to be repaired.

SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from COMECTA, S.A.

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Description

- Outlet filter Cod. 5900620
Outlet filter Cod. 5900621
Scale pressure gauge
Oil for vacuum pump GS-32, 1L

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of COMECTA, S.A.

Any manipulation of the apparatus by personnel not authorized by COMECTA, S.A. automatically cancels the guarantee.

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Para garantizar la seguridad del equipo los recambios deben adquirirse a COMECTA, S.A.